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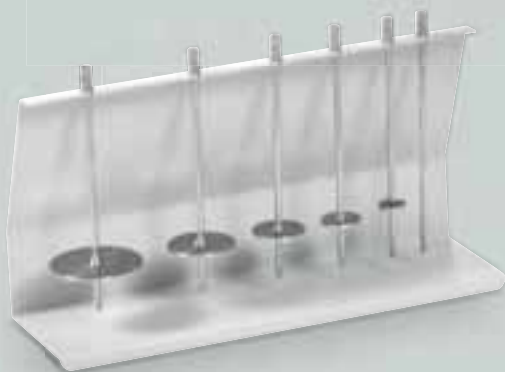
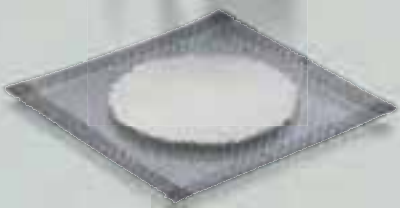
Intended use

- Raypa viscometers, models RP1 and RP2, are rotational viscometers for the quick determination of viscosity as specified in ISO 2555 and other ASTM norms.
- Instrument offers a precise digital torque measurement and viscosity measurement results which are 100% compatible to the Brookfield method.

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Features

- Interface RS232 for data transfer to PC or printer, temperature probe PT100, digital display for direct reading of various parameters, software Viscosoft Basic for easy data evaluation, warning of under- or over-range condition, complete spindle set and 24 months warranty.
- Additionally model RP2 offers bi-directional interface RS232 for RP2 control and data evaluation through a PC and the software Viscosoft Plus for the automated control of viscometer and the issuing of rheological studies.

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Compatibility

- ISO 2555 and ASTM define a viscometer in torque, speed and spindle design. Our Raypa viscometers comply with such specifications and are therefore 100% compatible to Brookfield method and able to perform measurements as per such norms.

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Additional information

- Our viscometers are supplied as complete equipment in a suitable carrying case. Standard accessories for each basic equipment: complete set of spindles with storage rack, interconnecting RS232 cable, spindle guard and instructions manual, are included in case.

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Displayed parameters

- Selected speed: rpm.
- Selected spindle: S.P.
- Dynamic viscosity: mPas o cP.
- % Scale (Brookfield): %.
- Sample temperature: °C o °F.
- Auto range to display viscosity limits: mPas o cP.
- Shear ratio S.R. (with special spindles. Only in V2 version): 1/s.
- Shear stress S.S. (with special spindles. Only in V2 version): N/m² o dyne/cm².

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Viscosity range

- Version L: 3 - 2000000 mPas in 76 ranges - 19 speeds with 4 spindles (for low to medium viscosity test fluids).
- Version R: 20 - 13000000 mPas in 114 ranges - 19 speeds with 6 spindles (for medium to higher viscosity test fluids).

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Supplied with

- Set of standard spindles, stand and carrying case.

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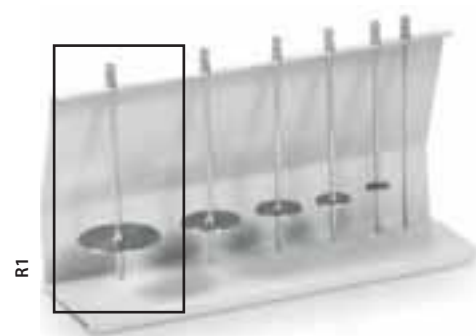
- 5 - 33300 mPas.

Spindle R1

- For an accurate measurement of low viscosity. Compatible with viscometers type R.

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Referencia: APM.

Features

- Accessory to allow accurate viscosity measurements of small sample volumes (from 8 to 13 ml).
- Works with all models.
- The sample chamber fits into a circulating chamber with controlled temperature. Using a suitable external thermostatic unit, sample can be thermostated between $-10\text{ }^{\circ}\text{C}$ and $+100\text{ }^{\circ}\text{C}$.
- It is supplied with a set of 3 or 4 spindles depending on version:
- Version L: TL5 - TL6 - TL7.
- Version R: TR8 - TR9 - TR10 - TR11.
- Viscosity range.
- Version L: from 1.5 to 200000 mPas/cP.
- Version R: from 25 to 3300000 mPas/cP.

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Features

- The "Adapter for Low Viscosity Materials" is necessary to obtain accurate measurements on low viscosity materials, both newtonian and no newtonian.
 - It can be used with V1 and V2.
 - It is supplied with a precision cylindrical spindle (LCP).
 - Reproducibility of viscosity measurements from 0.45 mPas.
 - Sample volume between 16-18 ml.
 - The "Adapter for Low Viscosity Materials" has a circulating chamber containing the sample tube. Through an external element (bath circulator), temperature in sample can be adjusted and controlled (temperature range between $-10\text{ }^{\circ}\text{C}$ and $+100\text{ }^{\circ}\text{C}$).
 - Viscosity range.
 - Version L: 0.3 to 2000 mPas/cP.
 - Version R: 3.2 to 21333 mPas/cP.
- Reference: ALV.**

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Features

- The adapter for helicoidal movement together with special T-shape spindles allows measurement of viscosity in substances as cream, gelatine, paste, wax, mastic, etc. Standard spindles cannot be used.
- The head automatic movement up and down avoids breaking of the sample structure to analyze, obtaining viscosity reliable readings.
- Height adjustable.
- Compatible with all models in our range of viscometers.
- It is supplied with: six suitable spindles with T-shape (PA, PB, PC, PD, PE, PF).
- Range of viscosity.
- Version L: from 156 to 3120000 mPas/cP.
- Version R: from 1660 to 33300000 mPas/cP.

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